

<b>160i or i160 Incubator</b>	Vios 160i and i160 Steri-Run Instructions	November 19, 2024

## Vios 160i and i160 Steri-Run Instructions

To run a Steri-Run decontamination routine on the Heracell Vios 160i, follow these steps:

1. **Prepare the Device:**
  - Ensure all samples and accessories (external probes, Hepa Filter, ect.) are removed from the workspace to avoid damage during the decontamination process. **Note: Unit Temp sensor stays in the unit**
  - Drain water from the reservoir and wipe up any residual water.
  - Make sure to remove the silicone plug from the inner chamber side and plug it into the outer side of the access port to avoid damage to the silicone.
2. **Start the Steri-Run:**
  - Press the Steri-Run key to display the Steri-Run menu.
  - Press the START key to activate the Steri-Run. The Steri-Run Progress dialog window will be displayed.
3. **Initial Steps:**
  - Open the glass door until the buzzer sounds (approximately 30 seconds).
  - Close the door after the buzzer sounds.
4. **Run the Steri-Run:**
  - The Steri-Run decontamination routine will start, and the device will heat up. The electromechanical door lock will engage at 65°C.
  - The display will show the current status, including temperature, start time, phase, and remaining run time.
5. **Completion:**
  - The routine will take approximately 12 hours to complete.
  - Once the routine is complete, open both device doors to air the workspace when the audible alarm sounds.
6. **Canceling Steri-Run (if needed):**
  - The Steri-Run decontamination routine can be interrupted at any time by pressing the Stop key. If the Stop key is pressed, the Steri-Run stop dialog box will be displayed for a safety scan.

Please refer to the user manual for any additional details or safety precautions.